

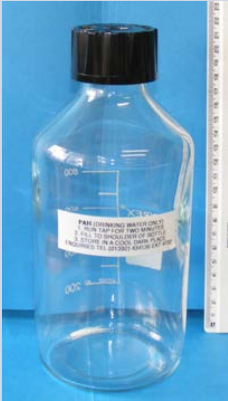
ORGANICS SAMPLE BOTTLES

BOTTLE USE	IMAGE	DESCRIPTION	SAMPLING COMMENTS	PRE - DOSING	MATRIX
Mass spec		2 L Screw cap glass	<ul style="list-style-type: none"> • Fill to neck • Dark and cold 	No dose	Clean waters
Neutral herbicides		100 ml Screw cap glass	<ul style="list-style-type: none"> • No rinse • Fill to shoulder • Cold 	1-2mg Sodium Sulphite	Clean waters

ORGANICS SAMPLE BOTTLES

BOTTLE USE	IMAGE	DESCRIPTION	SAMPLING COMMENTS	PRE - DOSING	MATRIX
Acid Herbicides		100 mL Screw cap glass	<ul style="list-style-type: none"> • No rinse • Fill to shoulder • Cold 	10-20mg Sodium Sulphite	Clean waters
Insecticides		1 L Screw cap glass	<ul style="list-style-type: none"> • No rinse • Fill to shoulder • Cold 	1.5 - 2.5g Sodium Thiosulphate	Clean waters

ORGANICS SAMPLE BOTTLES

BOTTLE USE	IMAGE	DESCRIPTION	SAMPLING COMMENTS	PRE - DOSING	MATRIX
Poly-aromatic-hydrocarbons		1 L Screw cap glass	<ul style="list-style-type: none"> • No rinse • Fill to shoulder • Cold 	10-20mg Sodium Sulphite	Clean waters
Organophosphorus pesticides and synthetic pyrethroids		500ml PET	<ul style="list-style-type: none"> • Fill to shoulder • Cold 	No dose	Clean waters
Hydrocarbons		1 L Screw cap glass	<ul style="list-style-type: none"> • Fill to shoulder • Dark & cold 	No dose	Clean waters

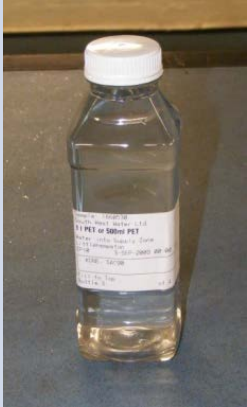

ORGANICS SAMPLE BOTTLES

BOTTLE USE	IMAGE	DESCRIPTION	SAMPLING COMMENTS	PRE - DOSING	MATRIX
Trihalomethanes		125ml SCREW CAP BROWN GLASS	<ul style="list-style-type: none"> • No rinse • Fill to top • Dark & cold 	150 - 250mg Sodium Thiosulphate	Clean waters
Geosmin, MIB & Metaldehyde		1 L SCREW CAP GLASS	<ul style="list-style-type: none"> • No rinse • Fill to shoulder • Cold 	No dose	Clean waters


ORGANICS SAMPLE BOTTLES

BOTTLE USE	IMAGE	DESCRIPTION	SAMPLING COMMENTS	PRE - DOSING	MATRIX
Oil type		500ml PET	<ul style="list-style-type: none"> • Dark & cold 	No dose	Clean waters
Total oil & grease in waste		1 L Screw cap glass	Screw cap glass	No dose	Waste waters
Volatile Acids		500ml Opaque plastic pot	<ul style="list-style-type: none"> • If sludge is very thin fill two pots 	No dose	Sludge

ORGANICS SAMPLE BOTTLES

BOTTLE USE	IMAGE	DESCRIPTION	SAMPLING COMMENTS	PRE - DOSING	MATRIX
Xylene & Toluene		250ml PET	<ul style="list-style-type: none"> • Fill to shoulder • Cold 	No dose	Various
Oils & fats		500ml 2 X Opaque plastic pot	<ul style="list-style-type: none"> • Please fill both pots 	<ul style="list-style-type: none"> • Fill to top • Cold 	Sludge

ORGANICS SAMPLE BOTTLES

BOTTLE USE	IMAGE	DESCRIPTION	SAMPLING COMMENTS	PRE - DOSING	MATRIX
NDMA		1 L PET	<ul style="list-style-type: none"> • Do not rinse • Fill to rim 	50mg Ascorbic Acid	Various

ORGANICS SAMPLE BOTTLES

BOTTLE USE	IMAGE	DESCRIPTION	SAMPLING COMMENTS	PRE - DOSING	MATRIX
Pentachlorophenol		50 ml Amber glass bottle wide neck	<ul style="list-style-type: none"> Cool & dark Return asap 	No dose	Various
Pentachlorophenol		500ml Amber glass bottle wide neck	<ul style="list-style-type: none"> Cool & dark Return asap 	No dose	Various
Permethrins		500ml Amber glass bottle wide neck	<ul style="list-style-type: none"> Cool & dark Return asap 	No dose	Various
Solvents		40ml Blue capped brown glass vial	<ul style="list-style-type: none"> Dark & cold 	Hydrochloric acid	Various
Organotins		1 L Screw cap glass	<ul style="list-style-type: none"> Fill to top Cold 	No dose	Various
Asulam		500ml Screw cap amber glass	<ul style="list-style-type: none"> Do not rinse Fill to rim 	0.5ml Sodium Thiosulphate solution	Various
Haloacetic acids		100ml Amber glass bottle	<ul style="list-style-type: none"> Fill to shoulder Cold 	0.1ml Ammonium Chloride solution	Various
Herbs		1 L Screw cap glass	<ul style="list-style-type: none"> Fill to top Cold 	No dose	Various